
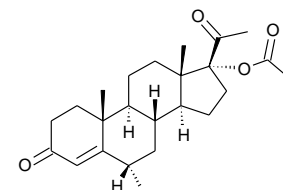


メドロキシプロゲステロン酢酸エステル(日本薬局方第十七改正記載条件)  
ーワイドポアC18カラムの活用ー


R191011A

## YMC-Triart Bio C18 (細孔径30 nm)

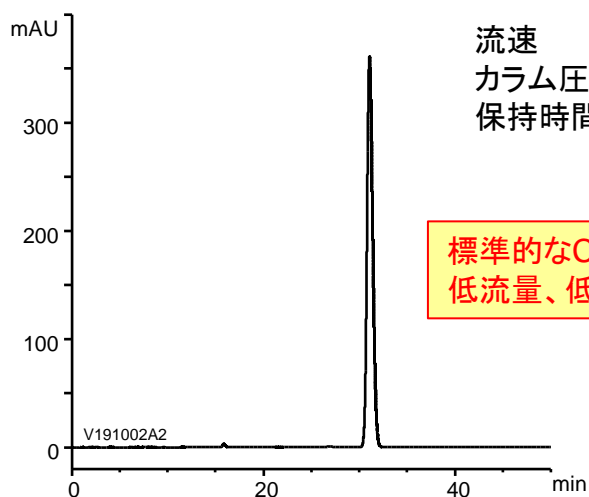

Medroxyprogesterone acetate

## YMC-Triart C18 (細孔径12 nm)

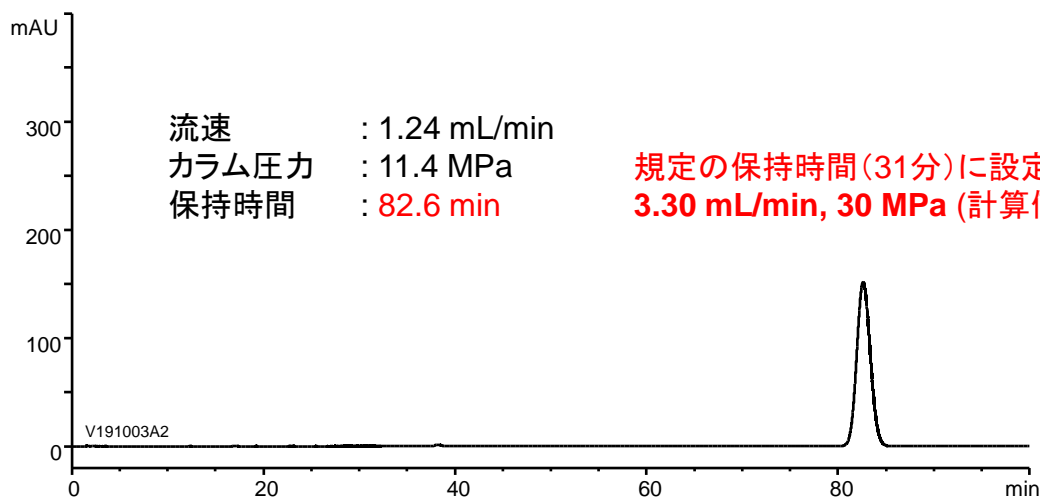

Column : 5  $\mu$ m, 250 X 4.6 mm I.D.  
Eluent : water/acetonitrile (3/2)  
Flow rate : 1.24 mL/min (adjust so that the retention time of medroxyprogesterone acetate is about 31 minutes)  
Temperature : 25°C  
Detection : UV at 254 nm  
Injection : 10  $\mu$ L  
(The Japanese Pharmacopoeia 17th; Assay)